nunchuck







Get sweet nunchuck skills


Nunchuck is the first and only platform that combines two key steps needed to make the perfectly-sized LNP – low-volume screening and scaled-up LNP formulation. Screen 8 total flow rates (TFRs) in just 10 minutes to find your sweet spot fast, then jump straight into making 0.250 mL to 1 L of your LNPs – all in the same day, on the same consumable. Dump the one-flow-rate-at-a-time, back-and-forth workflows for the ultimate one-and-done LNP maker.

0.250 mL-1 L formulations
Screen 8 TFRs in 10 min




The nun gets it done

Load your syringes filled with aqueous and organic solutions into Nunchuck, lock them into the Nun consumable, and you're ready to start making LNPs. The Nun's novel microfluidic mixing chambers can handle flow rates from 0.5 to 30 mL/min which means it can create a wide range of LNP sizes. Plus the nun can be used more than once – same day, same chemistry, same nun.

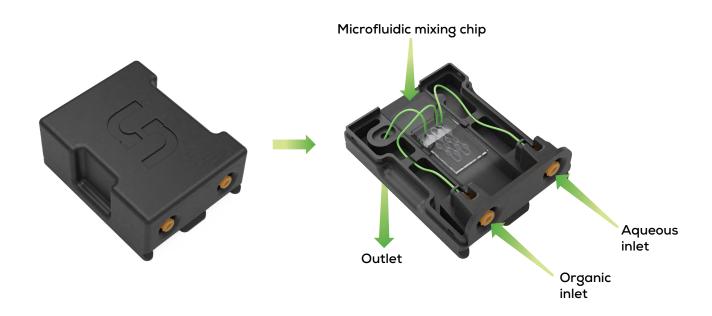


White belts welcome

Nunchuck is so easy to use a newbie can kick off a run in minutes. First, choose either a screening or prep run, then enter all the details about your formulation. The last step is to choose your flow rate ratio, total flow rates, and desired sample volume. You'll be an LNP black belt in 5 minutes flat.

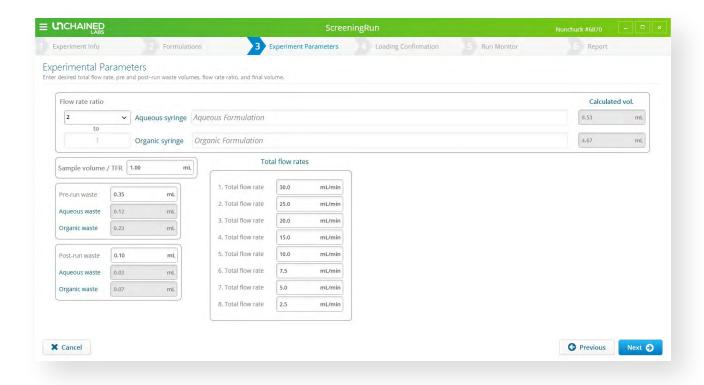


Hit your target

Depending on your target, you'll need to make different sizes of LNPs. Dialing in the total flow rate (TFR) is one of the most time-consuming steps for getting the size right. Nunchuck is outfitted with a fit-for-purpose screening run that nails down your TFR in 10 minutes. Screen up to 8 total flow rates with crazy repeatability so you can find your optimal conditions without breaking a sweat.

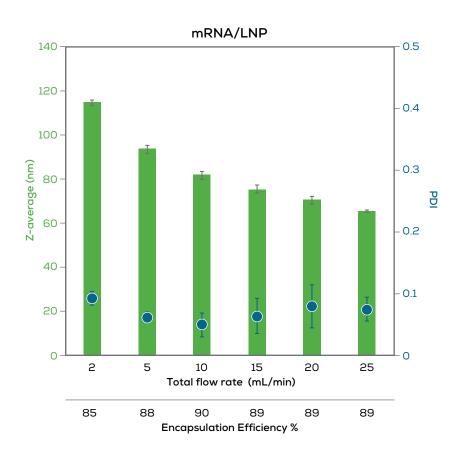

Level up

It doesn't matter if you need 0.250 mL to 1 L – Nunchuck gets you the same high-quality LNPs at any volume. Forget using one instrument to optimize your sample and another to make enough to characterize and develop your LNP formulations, Nunchuck always has your back.

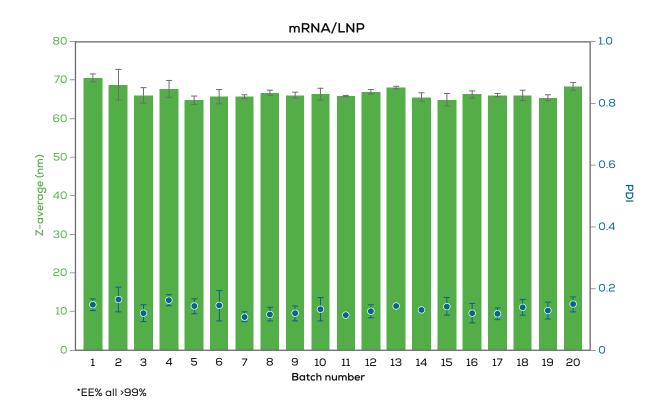

Specifications

Application	
Prep run volume range	0.250 mL-1 L
Flow rate ratios (aqueous to organic formulations)	1:1 to 10:1
Total flow rates (TFRs)	0.5-30 mL/min
Full LNP size range	50-200 nm*
Target polydispersity index	≤0.2*
Encapsulation Efficiency percent (EE%)	≥85%*
Instrument	
Environmental conditions	Temperature range: 18-28 °C Humidity: 40-60% relative humidity (non-condensing)
Physical	Weight: 26 kg Dimensions: 49 cm L x 40 cm D x 27 cm H
Physical Electrical	
,	Dimensions: 49 cm L x 40 cm D x 27 cm H

^{*} Sample and formulation dependent





Unchained Labs

6870 Koll Center Parkway Pleasanton, CA 94566 Phone: 1.925.587.9800 Toll-free: 1.800.815.6384 Email: info@unchainedlabs.com

© 2023 Unchained Labs. All rights reserved. The Unchained Labs logo, Nunchuck, and the Nunchuck logo are trademarks and/or registered trademarks of Unchained Labs. All other brands or product names are trademarks owned by their respective organizations.